

VORTEX-GENIE® 1 TOUCH MIXER

OPERATING INSTRUCTIONS

Models SI-0136 through SI-0186



70 Orville Drive, Bohemia, NY 11716 U.S.A.
(631) 567-4700 • Fax: (631) 567-5896 • Toll Free: 888-850-6208
Customer Service: custsvc@scientificindustries.com
www.scientificindustries.com

The latest Vortex-Genie® 1 Mixer is a powerful basic mixer that will easily obtain miscibility with the most difficult mixing situations. With proper care, the Vortex-Genie 1 will provide years of very reliable service and satisfaction. The Vortex-Genie 1 comes equipped with the standard Pop-off Cup attached. The Vortex-Genie 1 will also accept the 3-inch Platform.

APPLICATIONS

The standard Pop-off Cup will accommodate tubes of virtually any size or shape from micro-tubes to sizes that are 40mm in diameter and 200mm in length.



The 3-inch Platform accepts single or multiple test tubes, flasks, beakers, and other containers. The removable cover, made of thermoplastic rubber is highly impervious to chemical attack. It is recommended that spills are blotted quickly. Do not attempt to use the 3-inch Platform without its cover.



Catalog numbers:

OK-0500-902 3-inch Platform with Rubber Cover.
 OK-0500-901 3-inch Platform without Rubber cover.
 580-2013-00 Rubber cover for 3-inch Platform.

Vortex-Genie® is a registered trademark of Scientific Industries, Inc.
 U.S. Patent No. 4,781,487

OPERATING INSTRUCTIONS

1. Place the Vortex-Genie® 1 on a sturdy, level work surface.
2. Connect the power cord to a properly grounded three-pronged receptacle. 120V for the SI-0136 model; 230V for the SI-0146 model; 100V for the SI-0186 model.
3. The Vortex-Genie 1 is now ready for touch mixing. Simply touch a tube to the mixing cup, apply slight pressure, and the micro-switch will start the motor, thereby providing very vigorous mixing that will result in a vortex occurrence of the contents of the tube.
4. The 3-inch Platform is an accessory that may be ordered separately and used on the Vortex-Genie 1. The standard Pop-off Cup may be removed and replaced with the 3-inch Platform. **First unplug the Vortex-Genie 1 from its electrical power source.** Pop off the standard cup by inserting your fingers under the attachment flange and lift the Pop-off Cup with a prying motion. Align the flat surface of the Vortex-Genie 1 mixing shaft with the flat surface of the shaft receptacle on the under side of the 3-inch Platform. Press downward until the attachment snaps into place.
5. Reconnect the Vortex-Genie 1 to the properly grounded power source. The Vortex-Genie 1 is now ready for touch mixing using the 3-inch Platform.

⚠ CAUTION! Do not lift your Vortex-Genie 1 by its attachment.

Both the Pop-off Cup and the 3-inch Platform are removable. They will pop off easily if you lift them, and snap back on again just as easily by aligning the flats and pressing down.

⚠ CAUTION! Failure to follow operating instructions can compromise the user's safety.

Care and Handling

Your Vortex-Genie 1 Touch Mixer should be given the care normally required for any electrical appliance. Avoid wetting or unnecessary exposure to fumes. The finish can be washed with water (after unplugging) and soap or detergents, using a cloth or sponge. These and all replacement parts are available through most laboratory equipment distributors. Refer to "Parts Assembly List" and enclosed multi-fold brochure.

⚠ CAUTION! Unplug from power before cleaning. Do not immerse.

SPECIFICATIONS


The Vortex-Genie 1 Touch Mixer is classified as "Installation Category 2"
 Environmental: 0°C - 38°C (32°F-100°F), 95% Humidity max.

MODEL	POWER REQUIRED	AMPS
SI-0136	120V/60Hz.	0.65
SI-0146	230V/60Hz	0.5
SI-0186	100V/50-60Hz	1.0

VORTEX-GENIE® 1

PARTS ASSEMBLY LIST

To order parts for the Vortex-Genie 1:
 Contact your local distributor or visit
www.scientificindustries.com. Please specify
 Part No., quantity and electric voltage.

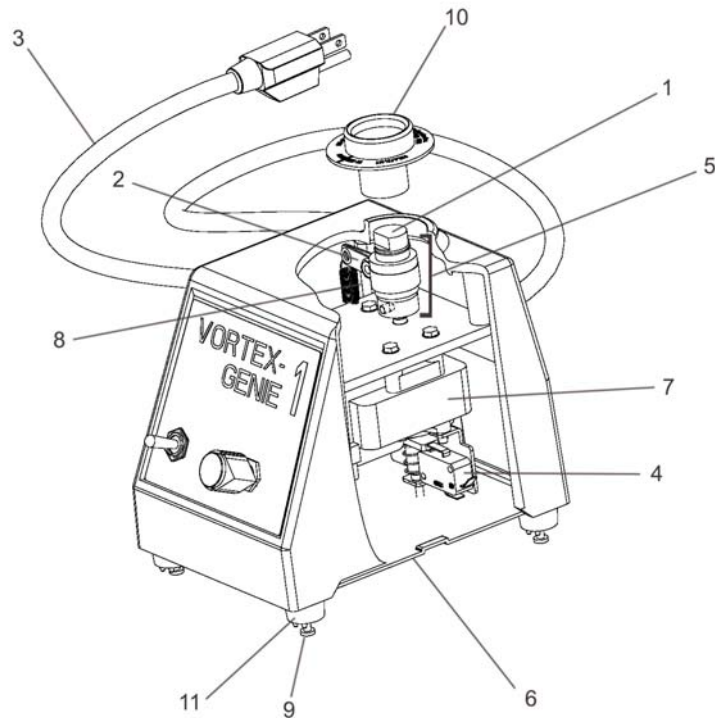

Indicator No.	Part No.	Description
1	OK-0236-902	Bearing Retainer Kit
2	OK-0236-903	Eccentric Clamp
3A	318-0510-02	120V Line Cord
3B	OK-0246-901	240V Line Cord, without Plug
3C	OK-0256-901	240V Line Cord, European Plug
3D	OK-0266-901	240V Line Cord, British Plug
3E	OK-0276-901	240V Line Cord, Swiss Plug
3F	OK-0286-901	100V Line Cord
4	ESP0005	Micro Switch
5	OK-0136-904	Eccentric with Clamp Assembly
6	0M-0236-210	Bottom Closure
7A	0A-0136-021	Motor 120V Assembly
7B	0A-0146-021	Motor 240V Assembly
7C	0A-0186-021	Motor 100V Assembly
8	0M-0136-201	Spring, Anti-Rotation
9	SCP0030	Screws, Bottom Closure (4)
10	146-3011-00	Pop-Off Cup
11	OK-0236-907	Feet with Screws (4)