

REPLACEMENT FEET WITH SCREWS

Part No. 0K-0236-900, Part No. 0K-0236-907,
Part No. 0K-0400-900, Part No. 0K-4000-900

For Models


SI-0236 through SI-0297	SI-D236 through SI-D297
SI-0136 through SI-0196	SI-T236 through SI-T296
SI-A236 through SI-A296	SI-0400 through SI-0406
SI-A536 through SI-A596	SI-4000 through SI-4006

INSTALLATION:

1. Unplug the mixer from any electrical outlet.
2. Remove any one of the bottom corner screws, leaving the remaining three in place.

NOTE: Do not remove all four screws to avoid exposing internal components.

3. Replace with the new screw and foot. Do not over tighten.
4. Repeat steps 2 and 3 for remaining three screws and feet.

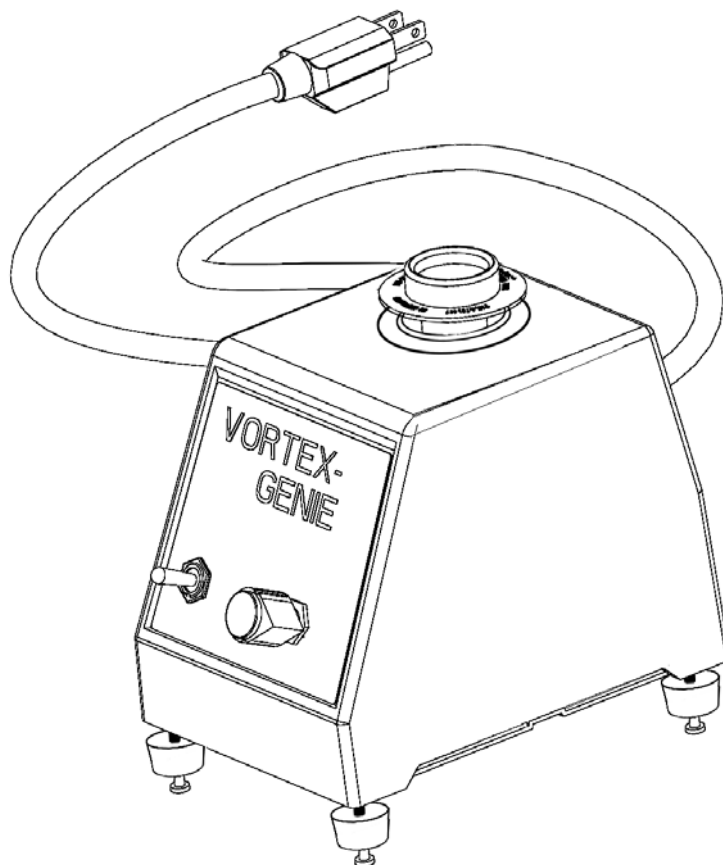

5. If your Vortex-Genie[®] 2 mixer has older style rubber strips for feet (one strip in the front and one in the back) these strips should be removed after the new feet are installed.



70 Orville Drive, Bohemia, NY 11716 U.S.A.
(631) 567-4700 • Fax: (631) 567-5896 • Toll Free: 888-850-6208
Customer Service: custsvc@scientificindustries.com
www.scientificindustries.com



© Scientific Industries Inc. 2011
LTP0017 Rev. E