

3-D ORBITAL SHAKER ATTACHMENT ACCESSORY FOR THE ENVIRO-GENIE®/INCUBATOR-GENIE™ Catalog No. SI-1250



1. Place 3-D Orbital Shaker Attachment inside the Enviro-Genie or Incubator-Genie chamber, aligning the hex shaft with the hex on the pulley of the 3-D Orbital Shaker Attachment. Make sure the two feet are in contact with the chamber base.
2. Use the Enviro-Genie/Incubator-Genie Rock Mode for controlling the speed of your 3-D Orbital Shaker Attachment. Speed range is 2-70RPM. Refer to the Enviro/Incubator-Genie manual for Rock Mode operation.
3. Place Non-Skid Mat on the 3-D Orbital Shaker platform. Do Not use attachment without Non-Skid Mat.
4. Place all loads centered on 3-D Orbital Shaker platform. Please note, platform capacity is 1 Kg (2.2 lb).



80 Orville Drive, Suite 102, Bohemia, NY 11716 U.S.A.
Phone: (631) 567-4700 • Fax: (631) 567-5896 Toll Free: 888-850-6208
Customer Service: custsvc@scientificindustries.com
www.scientificindustries.com