

ROTO-SHAKE GENIE®

STACKABLE NON-SKID TRAY

Operating Instructions

Model SI-1170



The Stackable Non-Skid Tray is an accessory for the ROTO-SHAKE GENIE®. The Stackable tray sits on top of the standard tray supplied with the ROTO-SHAKE GENIE® doubling the capacity of specimens being “rocked”.

OPERATING INSTRUCTIONS

1. Power “OFF” the ROTO-SHAKE GENIE. Place the Stackable Non-Skid Tray on top of the Non-Skid Tray supplied with the ROTO-SHAKE GENIE.
2. Power “ON” the ROTO-SHAKE GENIE with the SPEED set at “1”. Slowly ramp up to faster speed.



CAUTION! Do not start operation of the ROTO-SHAKE GENIE with the Stackable Non-Skid Tray at a high speed setting. Start at speed “1” and slowly amp up to faster speed.

CARE AND HANDLING

The ROTO-SHAKE GENIE Stackable Non-Skid Tray should be given the care normally required for any piece of laboratory equipment. The finish can be washed with a damp cloth and soap or mild detergents. Keep the item clean by immediately blotting any spills.



80 Orville Drive, Suite 102, Bohemia, NY 11716 U.S.A.
Phone: (631) 567-4700 • Fax: (631) 567-5896 Toll Free: 888-850-6208
Customer Service: custsvc@scientificindustries.com
www.scientificindustries.com