

ORBITAL SHAKER ATTACHMENT ACCESSORY FOR THE ROTO-SHAKE GENIE®


Installation and Operating Instructions


Model SI-1175



1. Remove magnetic platform from the Roto-Shake Genie® (see Roto-Shake Genie® manual).
2. Place the Orbital Shaker Attachment on the Roto-Shake Genie® base, aligning the hex shaft with the hex on the pulley of the Orbital Shaker Attachment. Make sure the two feet are in contact with the Roto-Shake Genie® base.
3. Use the Roto-Shake Genie® speed control knob for controlling the speed of your Orbital Shaker Attachment. Speed range is 15-100RPM.
4. Place all loads centered on Orbital Shaker platform. Please note, platform capacity is 2 Kgs (4.4 lbs).

 **CAUTION!** Remove from Roto-Shake Genie® before cleaning. Do not immerse.

 **CAUTION!** Do not use the Orbital Shaker Attachment in applications such as mixing of flammable materials or where transfer of mechanical energy to glass apparatus could lead to breakage..

 **CAUTION!** Failure to follow operating instructions can compromise the user's safety.

CARE AND HANDLING

The Orbital Shaker Attachment should be given the care normally required for any appliance. Avoid wetting or unnecessary exposure to fumes. The finish can be washed with a damp cloth and soap or mild detergents. Keep the unit clean by immediately blotting any spills.



80 Orville Drive, Suite 102, Bohemia, NY 11716 U.S.A.
Phone: (631) 567-4700 • Fax: (631) 567-5896 Toll Free: 888-850-6208
Customer Service: custsvc@scientificindustries.com
www.scientificindustries.com