

3-D ORBITAL SHAKER ATTACHMENT

ACCESSORY FOR THE ROTO-SHAKE GENIE®

Catalog No. SI-1150



1. Remove magnetic platform from the Roto-Shake Genie (see Roto-Shake Genie manual).
2. Place 3-D Orbital Shaker Attachment on the Roto-Shake Genie base, aligning the hex shaft with the hex on the pulley of the 3-D Orbital Shaker Attachment. Make sure the two feet are in contact with the Roto-Shake Genie base.
3. Use the Roto-Shake Genie speed control knob for controlling the speed of your 3-D Orbital Shaker Attachment. Speed range is 10-70RPM.
4. Place Non-Skid Mat on the 3-D Orbital Shaker platform. Do Not use attachment without Non-Skid Mat.
5. Place all loads centered on 3-D Orbital Shaker platform. Please note, platform capacity is 1 Kg (2.2 lb).



80 Orville Drive, Suite 102, Bohemia, NY 11716 U.S.A.
Phone: (631) 567-4700 • Toll Free: 888-850-6208
Customer Service: custsvc@scientificindustries.com
www.scientificindustries.com